


## 8 Channel Stopwatch

ordering code 10T6090

### Mode of operation:

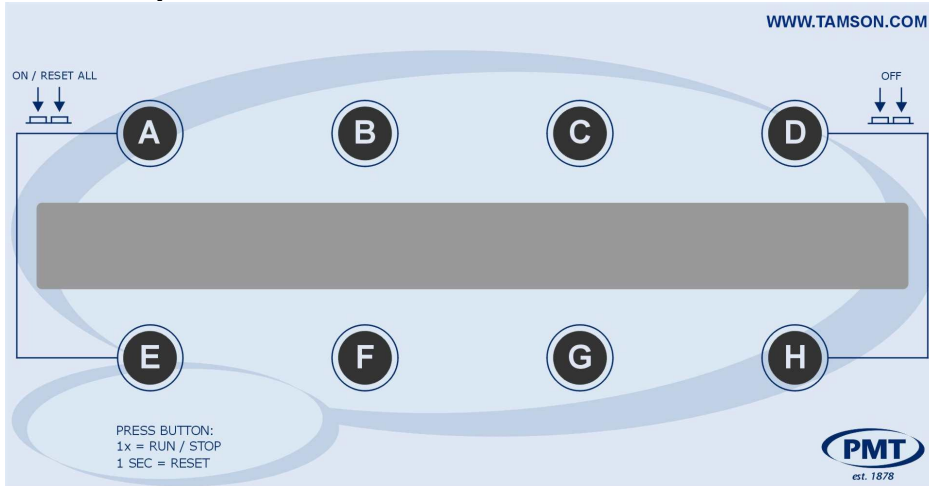

Eight individual stopwatches (A to H) each can be operated by it's individual switch:

- Press switch to start,
- Press switch to stop,
- Press switch to continue,
- Keep switch pressed for 2 seconds to reset stopwatch.

When power off, pressing one of the switches (A..H) will power-up the stopwatch.

Press both keys A+E to reset all stopwatches.

Press both keys D+H to switch the unit off.


When no timers are running the stopwatch automatically turns itself off, keeping current timing values in memory.

Replace battery when contrast dimms.

Battery life is minimum 6 months in normal mode of operation.

### Replacing battery:

Open the box at all 4 corners and lift top gently  
The 9V battery can easily be taken out

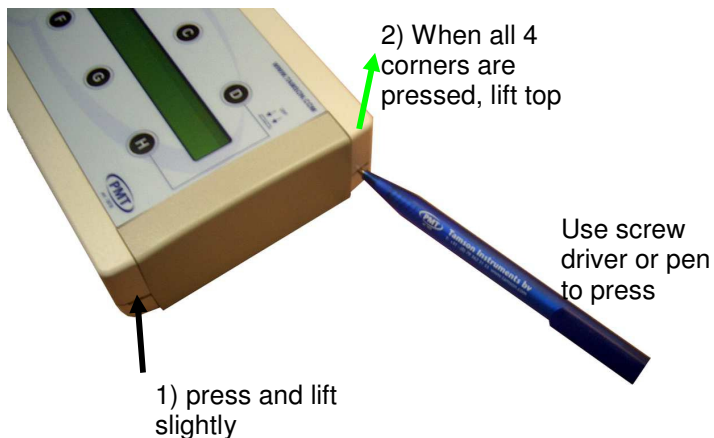

## Tamson Instruments bv